


Operation

1. Insert the sampler into the product. Ensure that the tip is inside the sampler body
2. At the required depth pull up the body to expose the tip. The sample will fall in around the tip
3. Push down body of the sampler to trap the sample
4. Withdraw sampler
5. Pull up body to empty sample

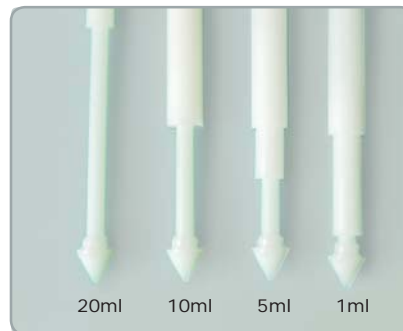

Available with a range of different tip sizes: 1ml, 5ml, 10ml and 20ml

| Stock No. | Pre-Sterile | Capacity | Length | Diameter | Material | Per Box |
|----------------|-------------|----------|--------|----------|----------|---------|
| 675 8060H501 | No | 1ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H501S | Yes | 1ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H505 | No | 5ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H505S | Yes | 5ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H510 | No | 10ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H510S | Yes | 10ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H520 | No | 20ml | 500mm | 21mm | HDPE | 20 |
| 675 8060H520S | Yes | 20ml | 500mm | 21mm | HDPE | 20 |
| | | | | | | |
| 675 8060H1001 | No | 1ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1001S | Yes | 1ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1005 | No | 5ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1005S | Yes | 5ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1010 | No | 10ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1010S | Yes | 10ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1020 | No | 20ml | 1000mm | 21mm | HDPE | 20 |
| 675 8060H1020S | Yes | 20ml | 1000mm | 21mm | HDPE | 20 |

QUICK QUOTE

ONLINE PRICING

Note: Other tip sizes are available between 1 and 20ml - subject to minimum order quantities
1500mm long samplers are available on request - subject to minimum quantities