

Disposable Slot Sampler

Cross Sectional Multilevel Sampler

The disposable Slot Sampler is ideal for taking a multilevel, cross sectional sample from powders and granules.

The Disposable Slot Sampler takes a single composite sample from the top, middle and bottom of your powder bed.


- Assembled & Packed in a Cleanroom
- Individually Bagged
- FDA & EU 10/2011 Conforming Materials
- BSE/TSE Free
- Full Batch Traceability
- ATEX Rated
- Available Sterile (gamma irradiated)



Operation

The Disposable Slot Sampler is quick and easy to use:

1. Push the sampler into the product
2. Rotate handle to open up slots
3. After leaving the sampler for several seconds to allow powder to enter, rotate handle to close the sampler.
4. Remove sampler
5. Slide the inner rod out of the outer tube
6. Deposit the sample into a suitable container

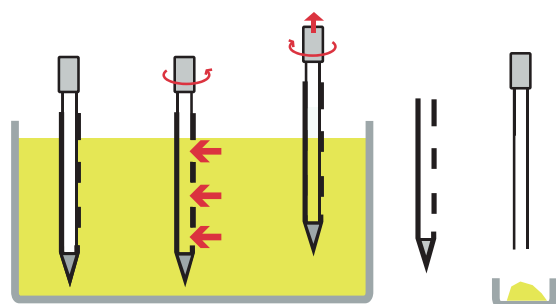

Stock No.	Pre-Sterile	Length	Vol	Diameter	Material	Per Box
675 8073H500	No	620mm	100ml	25mm	HDPE	20
675 8073H500S	Yes	620mm	100ml	25mm	HDPE	20
675 8073H1000	No	970mm	170ml	25mm	HDPE	20
675 8073H1000S	Yes	970mm	170ml	25mm	HDPE	20
6758073J1000	No	1000mm	500ml	40mm	HDPE	20
675 8073J1000S	Yes	1000mm	500ml	40mm	HDPE	20

[QUICK QUOTE](#)

[ONLINE PRICING](#)