

Micro Mate

Low Volume Powder Sample

The MicroMate is designed for low volume sampling of free flowing powders and granules.


The MicroMate can be easily adjusted so that it can take a range of sample sizes.

- Made from high quality 316 stainless steel
- Adjustable takes a range of sizes
- Simple to use





- 1. Insert the sampler into the product. Ensure that the inner rod extends beyond the outer tube.
- 2. At the required depth pull up the inner rod to draw the product into the sampler
- 3. Withdraw sampler.
- 4. Push down handle to release sample

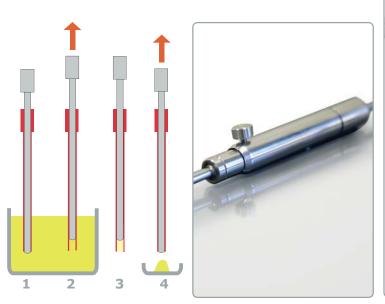





Stock No.	Sample Volume	Length	Material	Outer Diameter
675 1070A350	Variable - up to 0.1ml	350mm	316ss	5mm
675 1070A500	Variable - up to 0.1ml	500mm	316ss	5mm

QUICK QUOTE

ONLINE PRICING