


**Operation**

1. Insert the sampler into the product. Ensure that the tip is inside the sampler body
2. At the required depth pull up the body to expose the tip. The sample will fall in around the tip
3. Push down body of the sampler to trap the sample
4. Withdraw sampler
5. Pull up body to empty sample

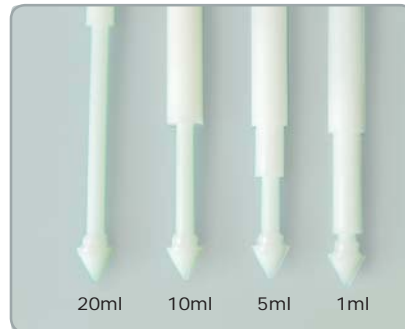

Available with a range of different tip sizes: 1ml, 5ml, 10ml and 20ml

**QUOTE REQUEST** ✓

**BUY ONLINE** ✓

Stock No.	Pre-Sterile	Capacity	Length	Diameter	Material	Per Box
675 8060H501	No	1ml	500mm	21mm	HDPE	20
675 8060H501S	Yes	1ml	500mm	21mm	HDPE	20
675 8060H505	No	5ml	500mm	21mm	HDPE	20
675 8060H505S	Yes	5ml	500mm	21mm	HDPE	20
675 8060H510	No	10ml	500mm	21mm	HDPE	20
675 8060H510S	Yes	10ml	500mm	21mm	HDPE	20
675 8060H520	No	20ml	500mm	21mm	HDPE	20
675 8060H520S	Yes	20ml	500mm	21mm	HDPE	20
675 8060H1001	No	1ml	1000mm	21mm	HDPE	20
675 8060H1001S	Yes	1ml	1000mm	21mm	HDPE	20
675 8060H1005	No	5ml	1000mm	21mm	HDPE	20
675 8060H1005S	Yes	5ml	1000mm	21mm	HDPE	20
675 8060H1010	No	10ml	1000mm	21mm	HDPE	20
675 8060H1010S	Yes	10ml	1000mm	21mm	HDPE	20
675 8060H1020	No	20ml	1000mm	21mm	HDPE	20
675 8060H1020S	Yes	20ml	1000mm	21mm	HDPE	20

Note: Other tip sizes are available between 1 and 20ml - subject to minimum order quantities  
1500mm long samplers are available on request - subject to minimum quantities