

Powder Thief


The PowderThief was originated over 20 years ago and is now a world leader in powder sampling. It is extensively used throughout the pharmaceutical, chemical and food industries around the world.

The PowderThief is suitable for sampling free flowing powders and granules. A range of tips is available so that a single sampler body can be used to accurately sample a range of different volumes.

- Made from high quality 316 stainless steel Wide range of sizes
- Simple to useO






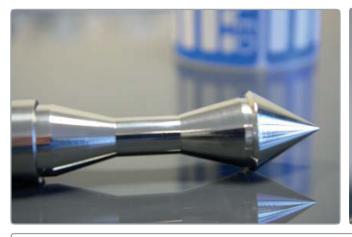



Operation

- 1. Insert the sampler into the product, ensure that the tip is inside the sampler body.
- At the required depth pull up the body to expose the tip. Powder will flow in around the tip.
- 3. Push down body of the sampler to trap the sample.
- 4. Withdraw sampler.
- Pull up body to release the sample.

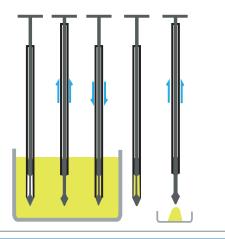





Sampler Bodies

QUOTE REQUEST \checkmark

•			
Stock No.	Length	Diameter	
675 1030A300	300mm	12.5mm	
675 1030A600	600mm	12.5mm	
675 1030A1000	1000mm	12.5mm	
675 1030A1500	1500mm	12.5mm	
675 1030A2000	2000mm	12.5mm	
675 1030B600	600mm	19mm	
675 1030B1000	1000mm	19mm	
675 1030B1500	1500mm	19mm	
675 1030B2000	2000mm	19mm	
675 1030C600	600mm	25mm	
675 1030C1000	1000mm	25mm	
675 1030C1500	1500mm	25mm	
675 1030D600	600mm	32mm	
675 1030D1000	1000mm	32mm	
675 1030D1500	1500mm	32mm	
675 1030D2000	2000mm	32mm	

Options Cleaning brushes See page 129
Storage cases See page 130
XRF Certification See page 139

Sampling Tips
A range of standard tip sizes is available for each body

BUY ONLINE

Stock No.	Vol.	Diameter		
675 1030AT025	0.25ml	12.5mm		
675 1030AT050	0.50ml	12.5mm		
675 1030AT075	0.75ml	12.5mm		
675 1030AT100	1.00ml	12.5mm		
/75 1000DT0		10		
675 1030BT2	2ml	19mm		
675 1030BT3	3ml	19mm		
675 1030BT5	5ml	19mm		
675 1030BT8	8ml	19mm		
675 1030BT10	10ml	19mm		
675 1030CT10	10ml	25mm		
675 1030CT20	20ml	25mm		
675 1030CT25	25ml	25mm		
675 1030DT20	20ml	32mm		
675 1030DT25	25ml	32mm		
675 1030DT50	50ml	32mm		
675 1030DT100	100ml	32mm		


... also available is a single use, disposable PowderThief

The Disposable PowderThief works in the same way as the stainless steel Powder Thief (above). Disposable samplers have the advantage that they can be used straight from the bag and never need to be cleaned, thereby saving you time and money.

See pages 12 & 13 for more information



Tel: (804) 318-3686

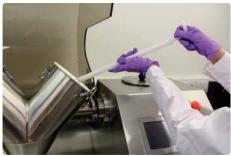



29