



**QAQC LAB**

www.qclabequipment.com


## Sleeve Sampler

Large Volume Sampling From Depth

Ideal for sampling large volumes at great depths. Various sized chambers can be attached to the same handle to allow different sample volumes to be taken.

- Single Point
- 316 stainless steel, bright finish
- Adjustable Length - use extension rods to make the sampler the required length

The heavy duty handle and extension rods have been designed to be extremely strong and robust to allow samples to be taken from depths up to 6m (after about 6m the sampler gets heavy).

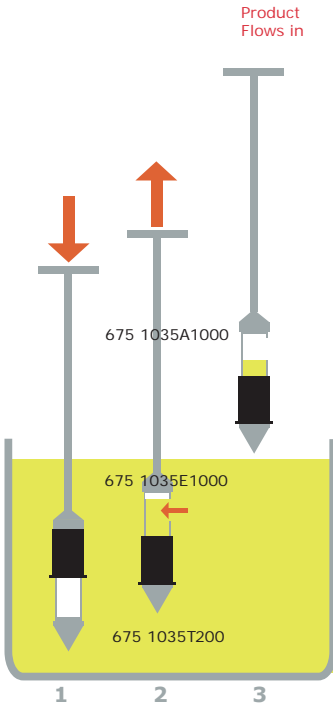

### Operation

1. Screw the Sample Chamber onto the handle and insert the sampler into the product.
2. At the required depth pull up the sampler slightly. This will force the sleeve down so that product can fall into the Sample Chamber.
3. Remove sampler.
4. To empty sampler either pour sample out of filling port or unscrew the tip.



[QUOTE REQUEST](#) ✓

[CLICK FOR PRICING](#) ✓

Handles & Extension Rods		
Stock No.	Description	Length
675 1035A1000	Heavy Duty Handle - 316ss	1000mm
675 1035A1500	Heavy Duty Handle - 316ss	1500mm
675 1035E1000	Heavy Duty Extension Rod - 316ss	1000mm
675 1035E1500	Heavy Duty Extension Rod - 316ss	1500mm

Sample Chambers		
Stock No.	Material	Chamber Volume
675 1035T100	316ss	100ml
675 1035T150	316ss	150ml
675 1035T200	316ss	200ml

Storage Cases	See page 129 for further information
XRF Certification	See page 139 for further information