


WHILE THE SAMPLE IS DRYING, TH RESULTS MAY BE VIEWED, AND LATER PRINTED AS A REPORT.

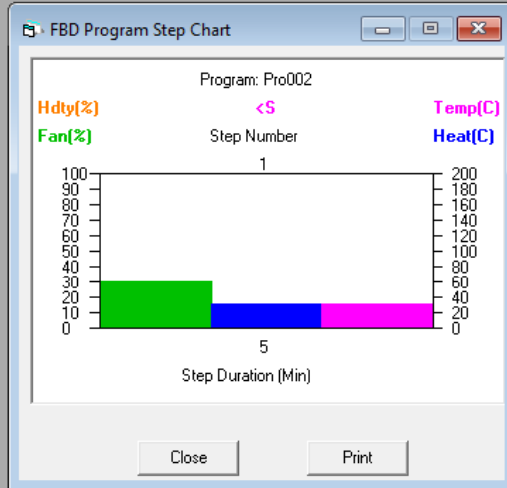

The "Control Program Manager" window shows settings for Program Pro002. It features a list of steps (1-16) on the left, with Step 1 selected and highlighted in green. The main area displays "Settings for Step 1 (Enabled)" under "Control parameters":

- Timer enabled: Timer duration (h:mm) 0:05
- Heater enabled: Heater set point 30.0°C
- Pause at end of step: Fan speed 30%

Below this is the "OR HUMIDITY" section with an "Auxiliary input test":

- Test enabled: Terminate: Step
- if Temperature is Less than 30.0°C

The "Pulser Unit" section shows:

- Enabled: Pulser valve closed 50%

At the bottom, there are buttons for "Execute", "Upload", "Save", and "Close". The "Steps" section includes "Show Chart", "Disable Step 1", "Insert after 1", and "Renumber".

DRYING VARIABLES MAY BE PROGRAMMED

OR PROGRAM